

## RGA OPTION 2 INSTALLATION INSTRUCTIONS

Prior to installing option 2 power supply, the RGA will have two identical skins. Only one of these skins needs to be removed to install the power supply. Looking at the rear of the RGA, the skin that is adjacent to the +24V DC external power connector is the one to be removed. Follow these easy steps.

1. Remove four screws on the skin to be removed. Save them for later, use when installing the option 2 power supply.
2. Remove skin and locate J601 header (towards rear of board).
3. Place connector from option 2 power supply on to J601 header. This connector can only be installed in one direction.
4. Install power supply such that the 120V AC power entry module faces the rear the RGA (away from the probe). The wires have extra length and must be tucked in so not to get pinched between the chassis and the power supply.
5. Install the four screws removed prior.

+24v dc  
external  
power  
connector



### STANFORD RESEARCH SYSTEMS

1290-D Reamwood Ave, Sunnyvale, CA 94089  
Phone (408) 744-9040, Fax (408) 744-9049  
e-mail: info@srsys.com, www.srsys.com