

Turn-On Procedure for the QMS Series Analyzers

1. With all four front-panel switches in the off position, turn on the power.
2. After the power has been turned on, turn on the Mechanical pump.
3. Now, turn on the Capillary Flow Valve.
4. When the pressure falls below 5 mbar, turn on the Turbo Pump. If a capillary is installed, it may take a little time to reach 5 mbar. Putting your finger over the end of the capillary will speed the process up.
5. When the Turbo Pump has reached full speed, turn on the Sample Inlet Valve.